

Instructions for use

Rotor A-2-VC

English (EN)

This document complements the operating manual for the Concentrator/Vacufuge® and does not replace it. Therefore, please also read the operating manual for the Concentrator before starting up the rotors for the first time. You can find the current version on the Internet at www.eppendorf.com.

1 Safety notice




NOTICE! If handled incorrectly, the rotor can fall over.

The rotor can fall over if the buckets are used as a handle.

- ▶ Remove the buckets before inserting and/or removing a swing-bucket rotor.
- ▶ Always use both hands to carry the rotor cross.


2 Inserting skirted plates

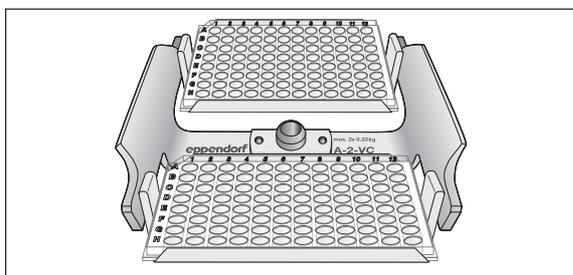

- ▶ Insert the microplates and PCR plates (skirted) into the buckets. Do not allow the plates to protrude over the edge of the bucket.

3 Inserting tubes, PCR strips and plates


Insert tubes, PCR strips and unskirted plates using the work plate (order no. 0030 124.235).

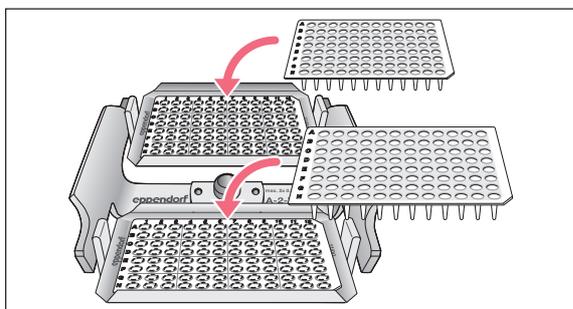

- ▶ Load the work plate.
- ▶ Fit the work plate into the cutout on the lower edge of the bucket using the guide.

4 Maximum filling

During centrifugation, the menisci in the tubes along the edges of the plates are at an angle.

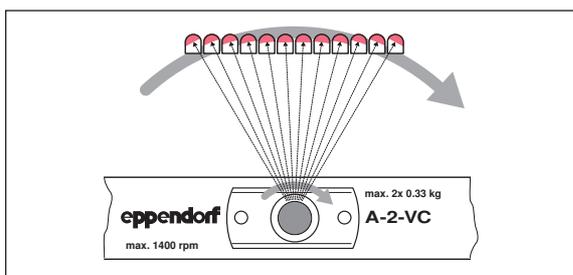

- ▶ Fill the wells of the plates with $\frac{2}{3}$ of the max. filling volume.

Your local distributor: www.eppendorf.com/contact

Eppendorf AG · 22331 Hamburg · Germany

eppendorf@eppendorf.com · www.eppendorf.com