

## Safety notice

Rotor S-4xUniversal (Centrifuge 5910 R)

English (EN)



This document is a supplement to the manual of the Centrifuge 5910 R and does not replace it. Therefore, please also read the manual of the centrifuge before starting up the rotors and accessories for the first time. The current version of the manual for all available languages can be found on the Internet under [www.eppendorf.com/manuals](http://www.eppendorf.com/manuals).



### **NOTICE! Material damage due to incorrect equipping of the swing-bucket rotor.**

Incomplete equipping of the swing-bucket rotor or an uneven load will reduce the life span of the rotor and the corresponding buckets considerably.

- ▶ Always load all positions of a swing-bucket rotor with buckets.
- ▶ Load opposite buckets with the same weight (adapter, tubes, or plates and content).

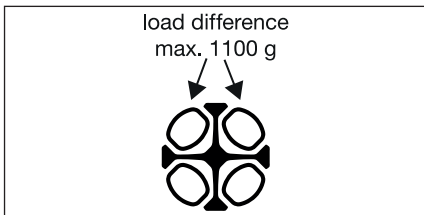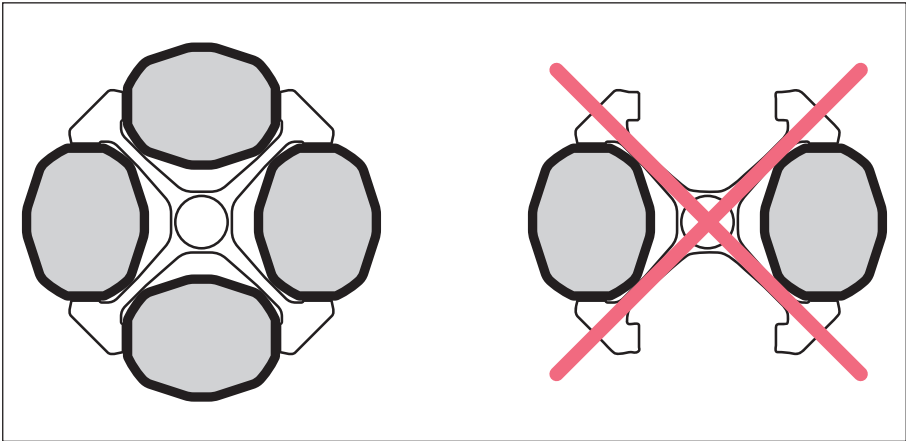

- ▶ Load adjoining buckets with a maximum of 1100 g difference in weight.

Your local distributor: [www.eppendorf.com/contact](http://www.eppendorf.com/contact)

Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany

[eppendorf@eppendorf.com](mailto:eppendorf@eppendorf.com) · [www.eppendorf.com](http://www.eppendorf.com)