
eppendorf

NEW BRUNSWICKTM SHAKER OPERATING KEY M1190-9600 Revision C

- 1. RPM: Lights when display shows speed, i.e., actual speed or setpoint in SET mode. Flashes when 5 RPM above or below setpoint.
- 2. HRS: Lights when display shows time, i.e., actual time remaining, or setpoint in SET mode, or total running time.
- 3. °C: Lights when display shows temperature, i.e., actual chamber or bath temperature, or setpoint in SET mode. Flashes when 1 °C above or below setpoint.
- 4. *: Innova 3100 Flashes when water level is below setpoint.
- 5. Determines the Function display: RPM, HRS, °C

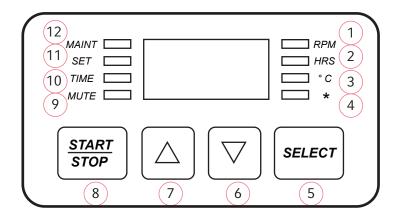

- 6. Decreases setpoint for agitation speed (RPM), time (HRS) or temperature (°C).
- 7. Increases setpoint for agitation speed (RPM), time (HRS) or temperature (°C).
- 8. Starts or stops shaking.
- 9. MUTE: Lights when audible alarm is deactivated.
- 10. TIME: Lights during a timed run. Flashes at end of run.
- 11. SET: For establishment of setpoints, lights when using \triangle or ∇ key
- 12. MAINT: Every 10,000 hours of operation, lights to signal need for preventative maintenance

Audible alarm:

- 1. Press SELECT to acknowledge. Stops sounding, but remains enabled.
- 2. To deactivate, SELECT to HRS, then press ▲and ▼ SIMULTANEOUSLY, then press START/STOP. MUTE will light.
- 3. To reactivate, SELECT to HRS, then press ▲ and ▼ SIMULTANEOUSLY, then press START/STOP. MUTE will go off.

To set and start a timed run:

- 1. Press SELECT until HRS lights.
- 2. Press ▲or▼to set time (between 0.1 HRS and 99.9 HRS).
- 3. While SET is lit, press START/STOP. TIME will light and Display will show time remaining.

To cancel:

- 1. Press SELECT to HRS.
- 2. Press ▲or ▼, then START/STOP. Time light turns off. Display reads "Off".

To adjust running time:

While TIME is lit, press ▲ or ▼ to set a new time.

Total running time:

- 1. Press SELECT to HRS.
- 2. Press ▲ and ▼ SIMULTANEOUSLY.
- 3. Press START/STOP. Display indicates hundreds of hours.

Flashing display indicates power interruption. Press any key to deactivate.